## NRPI radon chamber



### NRPI thoron/radon chamber



### I. Radon chamber

# **Details of calibration facility:**

**Typical radon** concentration:  $(100 - 8000) \text{ Bq/m}^3$ 

**Volume:** 45 m<sup>3</sup>, type: dynamic -Walk in, walls similar to dwellings

Continuously monitored radon gas, EEC, equilibrium factor,  $f_p$ , temperature, relative and controlled humidity, air exchange rate, aerosol concentration and size

**parameters:** distribution (5 - 1100) nm GRIMM (SMPS + C)

Available for solid (carnauba wax) polydisperse aerosols or liquid monodisperse

monitoring: (NaCl) aerosols.

**Uncertainties (95% CI):** radon gas < 10 %, EEC < 20 %.

# Typical adjustable and stable parameters in NRPI atmosphere:

Quantity	Typical used values
temperature	(8-45) °C
rel. humidity	(20 - 95) %
mean solid aerosol aerodynamic diameters	(100 – 200 ) nm
monodisperse liquid aerosol aerodynamic diameters	(0,5 - 3) μm
Air exchange rate	$(0,1-2) h^{-1}$
Equilibrium factor F	(1-90) %
$f_{ m p}$	(1-60) %

#### II. Thoron/radon chamber

## **Details of calibration facility:**

**Thoron concentration:**  $100 - 100\ 000\ \text{Bq/m}^3$ 

**Radon concentration:**  $0.1 \text{ Bq/m}^3 - 10 \text{ MBq/m}^3$ 

**Ratio** (K) of radon/thoron: arbitrary ratios in wide range: K < 1 or K > 1

**Volume:** 150 dm<sup>3</sup>, stainless steel vessel

**Continuously monitored** 

and stable parameters: radon gas, thoron gas, relative humidity, room temperature

**Uncertainties (95 CI):** radon gas < 10 %, thoron gas < 15%.

### Typical adjustable and stable parameters in NRPI atmosphere

Quantity	Typical used values
Temperature	(room temperature)
Rel. humidity	(5-60) %
$f_p$	> 50%

All of the NRPI reference instruments are independently compared with a primary instruments of the renowned Labs. as the PTB Braunschweig (Germany), BfS Berlin (Germany) and Czech Authorized Metrological Centre Kamenna.

Description of the QA Programme for radon gas and short lived radon progenies can be found here - doi:10.1093/rpd/ncn113 in the Radiation Protection Dosimetry, Vol. 130, No. 1, pp. 43-47, 2008.

#### **REFERENCES:**

#### 2015

2014 ICHLNRRA Intercomparison of Radon/Thoron Gas and Radon Short-Lived Decay Products Measuring Instruments in the NRPI Prague, doi:10.1093/rpd/ncv311

#### 2014

International Intercomparison of Measuring Instruments for Radon/Thoron Gas and Radon Short-Lived Daughter Products in the NRPI Prague, doi:10.1093/rpd/ncu079

#### 2011

International Intercomparison of Measuring Instruments for Radon Gas and Radon Short-Lived Daughter Products in the NRPI Prague, doi:10.1093/rpd/ncu079